

Environmental Swab

ETS Environmental Monitoring Swab



REST™ ETS Environmental KIT (Surface Inspection Kit)

- 1 By using assembled swab with cap and exclusive tube, it is easy and fast.
- 2 This product is for surface inspection of many kinds of instruments, device, human body and etc.
- 3 Various solution and volume of saline, BPW, Lethen Broth, NRS, Bufferfield solution, Dey/Engley Neutralising Broth.
- 4 Multiple swab applicators based on your target specimens.
- 5 Used for environmental monitoring of food, cosmetic, pharmaceutical and other facility.



REST™ ETS Swab How to use



- [Fig.1] Open the cap and push the swab head to the surface of inside wall to remove extra solution in the swab head.
- [Fig.2] Collect specimen by using 10x10 cm or 4x5 cm sample template
Sample operator needs training to obtain specimen to test.
- [Fig.3,4] After appropriate sample collection, put the swab into the tube and close the cap tight.
- [Fig.5] Transfer the tube containing specimen to laboratory for analysis as soon as possible. Transport sample tube in 24 hrs at room temperature and store nutrient Solution (Peptone, Lethen) at 4°C.

REST™ ETS Swab [Swab Applicator Type]



1. Standard Nylon Swab : General, Standard Swab Kit
2. Big head Nylon Swab : Larger head than standard (bigger surface)
3. Standard Flocked Swab : General, Standard Flocked Swab Kit
4. Mini tip Flocked Swab : Mini tip size and flexible stick



Environmental Swab

ETS Environmental Monitoring Swab




REST™ ETS Environmental Swab Media & Swab Applicator Type

1. Media Selection : Saline 0.85%, Buffered Peptone Water, Lethen Broth, Butterfield's Solution, Neutralising Rinse Solution, Dey/Engley Neutralising Broth
2. Media Volume : 1 ml, 5 ml, 10 ml and customer ordered
3. Swab Applicator : Standard Nylon Swab, Big head Nylon Swab, Standard Flocked Swab, Mini tip Flocked Swab

Ordering Information

Packaging : 50ea/Case

Component	Volume	Standard Nylon Swab Cat.No	Big head Nylon Swab Cat.No	Stadard Flocked Swab Cat.No	Mini tip Flocked Swab Cat.No
Saline 0.85%	5 ml	ETS-01	ETS-02	ETS-03	ETS-04
Saline 0.85%	10 ml	ETS-1	ETS-2	ETS-3	ETS-4
BPW (Buffered Peptone Water)	5 ml	ETS-05	ETS-06	ETS-07	ETS-08
BPW (Buffered Peptone Water)	10 ml	ETS-5	ETS-6	ETS-7	ETS-8
Lethen Broth	1 ml	ETS-10	ETS-13	ETS-16	ETS-19
Lethen Broth	5 ml	ETS-11	ETS-14	ETS-17	ETS-20
Lethen Broth	10 ml	ETS-12	ETS-15	ETS-18	ETS-21
Butterfield's Solution	5 ml	ETS-30	ETS-32	ETS-34	ETS-36
Butterfield's Solution	10 ml	ETS-31	ETS-33	ETS-35	ETS-37
NRS (Neutralising Rinse Solution)	1 ml	ETS-40	ETS-43	ETS-46	ETS-49
NRS (Neutralising Rinse Solution)	5 ml	ETS-41	ETS-44	ETS-47	ETS-50
NRS (Neutralising Rinse Solution)	10 ml	ETS-42	ETS-45	ETS-48	ETS-51
Dey/Engley Neutralising Broth	5 ml	ETS-60	ETS-62	ETS-64	ETS-66
Dey/Engley Neutralising Broth	10 ml	ETS-61	ETS-63	ETS-65	ETS-67

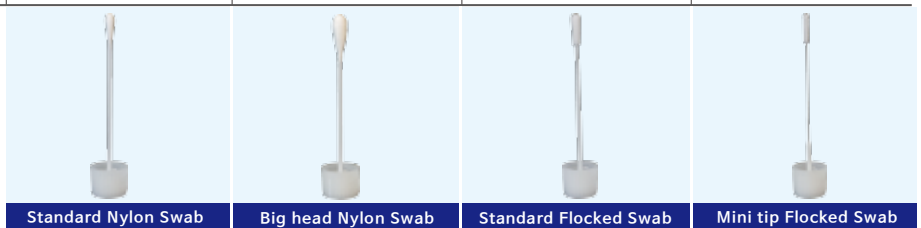

REST™ Diluent (9 ml Sterile Diluent(Saline and Phosphate Buffer))

Component	Cat.No	Packaging	Volume
Saline 0.85% Diluent	ETS-100	50 ea/Case	9 ml
Butterfield's Phosphate Buffered Diluent	ETS-101	50 ea/Case	9 ml



* Gamma-Irradiated Diluent with Certificate

